

### Summary and explanation of the kit:

OMNImet™-GUT provides the materials and instructions for the collection and stabilization of metabolites from human fecal samples.









### Warnings and precautions:

- Handle with care.
- Keep out of reach of children.
- Keep away from heat sources, sparks and open flames.
- Keep stabilizing liquid tube tightly closed when not in use.
- FOR EXTERNAL USE ONLY.
- Do NOT remove the yellow tube top from the tube.
- Do NOT spill the stabilizing liquid in the tube.
- Wash with water if liquid comes in contact with eyes or skin. Can cause serious eye irritation. Do NOT ingest.
- See MSDS at [www.dnagenotek.com](http://www.dnagenotek.com).
- If collecting a liquid stool sample, see separately provided user instructions.
- Small items may pose a choking hazard.

**Storage:** Room temperature storage pre- and post-collection 15°C to 25°C (60°F to 78°F).

Ship in accordance to applicable regulations covering transport of biological specimens.

### Label legend:

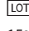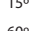-  Flammable
-  Warning
-  Collect sample by (Use by)
-  Catalog number
-  Manufacturer
-  Caution, consult instructions for use
-  Lot number
-  15°C / 25°C Storage instructions  
60°F / 78°F


### For Research Use Only

Not for use in diagnostic procedures

## USER INSTRUCTIONS

Read all instructions prior to collection.

### Procedure:

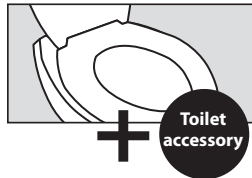

- 1 IMPORTANT PREPARATIONS:**
- Empty your bladder before beginning the collection.
  - Collect fecal sample free of urine or toilet water.
  - Utilize any toilet accessory which may be supplied with the collection device.
  - Toilet paper or tissues may be required.



- 2** While holding the yellow tube top, unscrew ONLY the purple cap and set aside for later use.
- Handle the stabilization liquid with extra care.**




**IMPORTANT:**  
Do NOT remove the yellow tube top.  
Do NOT spill the stabilizing liquid in the tube.  
Tube is not designed to stand upright, do NOT spill the stabilizing liquid in the tube.



- 3** Use the spatula to collect a small amount of fecal sample.



- 4** Transfer the fecal sample into the top portion of the yellow tube top.
- Repeat until the sample fills the yellow tube top.**

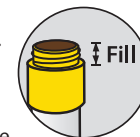

**IMPORTANT:** Do NOT push sample into the tube.




- 5** Scrape horizontally across the tube top to level the sample and remove any excess.
- Wipe exterior of tube and top with toilet paper or tissue as needed.



- 6** Pick up the purple cap and screw onto the yellow tube top until tightly closed.



- 7** For a **minimum of 30 seconds**, shake the sealed tube as hard and fast as possible in a back and forth motion.
- The fecal sample will be mixed with the stabilizing liquid in the tube; not all particles will dissolve.

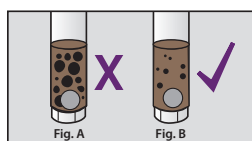

- 8** **IMPORTANT:** Continue shaking if large particles remain as shown in Figure A.



- 9** Place spatula in original packaging or wrap in toilet paper and discard in garbage.
- IMPORTANT:** Send the sample for processing following the delivery instructions supplied separately by the kit provider.