EN English – Instructions for sample collection.....	2
FR Français – Instructions de prélèvement de l'échantillon	3
ES Español – Instrucciones para la recogida de la muestra	4
DE Deutsch – Anweisungen zur Probenahme	5
NB Norsk – Instruksjoner for prøveinnsamling.....	6
SV Svenska – Anvisningar för provinsamling.....	7
DA Dansk – Instruktioner til indsamling af prøve.....	8
NL Nederlands – Instructies voor monstername	9
CS Český – Pokyny k odběru vzorku	10
IT Italiano – Istruzioni per la raccolta dei campioni.....	11
PL Polski – Instrukcje dotyczące pobierania próbeki	12
PT Português – Instruções para a recolha de amostras.....	13
BZ Português – Brasil – Instruções para coleta de amostras.....	14
AR العربية – تعليمات جمع العينات	15
ZH 中文 – 样本采集说明	16
KO 한글 – 샘플 채취 지침.....	17
JA 日本語 – サンプル採取に関する指示書.....	18
BG Български – Инструкции за вземане на проба	19
HR Hrvatski – Upute za prikupljanje uzoraka	20
ET Eesti keel – Juhised proovivõtmiseks.....	21
FI Suomi – Ohjeet näytteen ottamiseen.....	22
EL Ελληνικά – Οδηγίες συλλογής δείγματος.....	23
HU Magyar – Mintagyűjtési utasítások	24
IE Gaeilge – Treoracha maidir le bailiú an sampla	25
LV Latviešu valoda – Norādījumi parauga paņemšanai	26
LT Lietuvių k. – Mėginio paėmimo instrukcijos	27
MT Malti – Istruzzjonijiet għall-ġbir ta' kampjun	28
RO Română – Instrucțiuni pentru colectarea probelor.....	29
RU Русский – Инструкции для сбора образца	30
SK Slovensky – Pokyny na odber vzorky	31
SL Slovenščina – Navodila za jemanje vzorca:.....	32

Collection precautions:

Ensure the sponge tip does NOT come into contact with any surface prior to collection.

Donor should NOT eat, drink, smoke, or chew gum for 30 minutes before collecting saliva sample.

Intended use: For the collection and stabilization of RNA from human saliva samples.

Contents: 1 kit containing 1 mL stabilizing liquid.

Warning and precautions: Choking hazard. Caution should be used when inserting sponge into the mouth.

Wash with water if stabilizing liquid comes in contact with eyes or skin. Do NOT ingest. See SDS at www.dnagenotek.com.

Report any serious incident to DNA Genotek and the competent authority in your country.

Storage: 15°C / 25°C (59°F / 77°F)

Summary and explanation of the kit:

ORAcollection™ RNA is a self-collection kit that provides the materials and instructions for collecting human saliva samples.

Label legend:

RUO

Research Use Only

REF

Catalog number

Manufacturer

Do not reuse

Consult instructions for use

Caution

Use by

LOT

Lot number


QTY

Quantity of units per package


15°C / 25°C
59°F / 77°F

Storage instructions


MADE IN US
FROM FOREIGN
AND DOMESTIC
MATERIALS

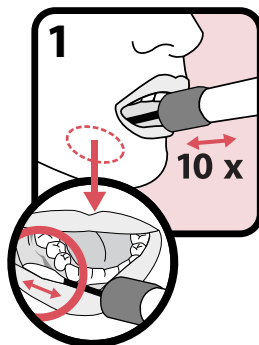Made in US from foreign
and domestic materials

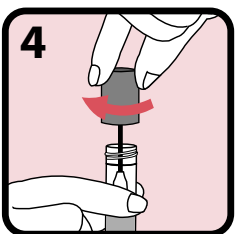
Open package and remove collector without touching sponge tip. Place sponge as far back in the mouth as comfortable and rub along the lower gums (see close-up image) in a back and forth motion. Gently rub the gums 10 times. If possible, avoid rubbing the teeth.

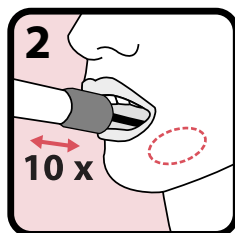
Gently repeat rubbing motion on the opposite side of the mouth along the lower gums for an additional 10 times.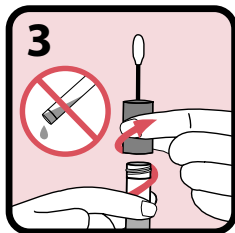

Hold the tube upright to prevent the stabilizing liquid inside the tube from spilling. Unscrew the cap from the collection tube without touching the sponge.

Turn the cap upside down, insert the sponge into the tube and close cap tightly.

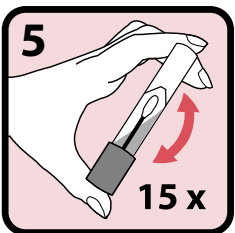Invert the capped tube and shake vigorously 15 times.

RUO

Research use only

DNAGENOTEK™

Superior samples
Proven performance

DNA Genotek Inc.
3000 - 500 Palladium Drive
Ottawa, ON, Canada K2V 1C2

Toll-free (North America): 1.866.813.6354
Tel.: +1.613.723.5757 • Fax: +1.613.723.5057
info@dnagenotek.com
www.dnagenotek.com

ORAcollection and DNA Genotek are trademarks of DNA Genotek Inc., registered under various jurisdictions. Some DNA Genotek products may not be available in all geographic regions; contact your sales representative for details. All DNA Genotek protocols, white papers and application notes are available in the support section of our website at www.dnagenotek.com.

Patent (www.dnagenotek.com/legal/notice)
© 2024 DNA Genotek Inc., a subsidiary of OraSure Technologies, Inc., all rights reserved.
PD-PR-1048 v2/2024-12

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ORAcollection in DNA Genotek sta blagovni znamki DNA Genotek Inc.
© 2024 DNA Genotek Inc., podružnica OraSure Technologies, Inc., vse pravice pridržane.